



TIPS: Filling the Spin Disk


1. Fill disk with syringe upright using slow, even pressure while depressing the syringe plunger. The lower level chamber will begin to fill with water and then the upper level chamber will fill with water. Fill the upper level chamber until the water level is between the two fill lines.

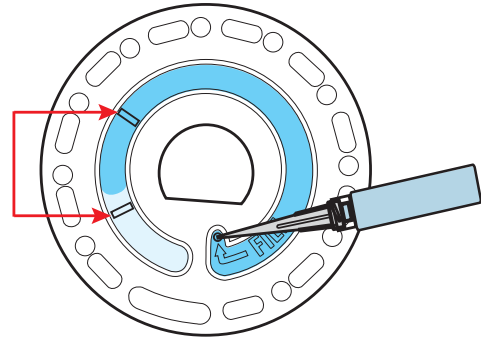

2. If the plunger is depressed too quickly an air bubble may develop half way through the filling procedure as shown in the diagram.

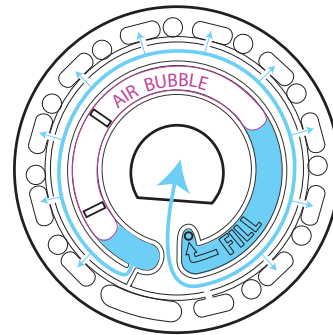

3. If an air bubble occurs, pick up the disk from the sides, tilt to the right side and rotate clockwise slowly until you displace the bubble.

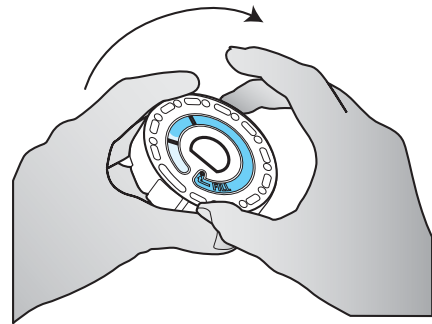

4. If the disk then shows a proper fill, between the fill lines, go ahead and use. If it is under-filled, add more of the water sample until it's between the fill lines.

